PHARMA
DIVISION

Pharmaceutical Industry
Equipment and Services

PATGROUP

Comasa

Wynka

Ecovida

Copharm

We are a multidisciplinary group of enterprises, specialized in equipment manufacture and productive projects analysis, advice and implementation. We offer comprehensive solutions to our clients, working with them in the entire process and promoting synergy through the joint work of the companies that make us up.

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PATGROUP

We design and build high-precision and quality machinery for the compression of powders and granules, as well as complementary process equipment.

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We design and build machinery of high precision and quality for powders compression, as well as complementary equipment of process.

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We provide services and integrate machines and process, according to customer and product needs. We close the circle between EQUIPMENT DESIGN, PRODUCTION PROCESS and PRODUCTION EFFICIENCY

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We are an engineering and construction company that is responsible for providing support regarding the location of machinery in the areas and the adaptation under current sanitary regulations. We also offer assembly and piping services completing a safe and efficient installation process.

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Turnkey Proposal

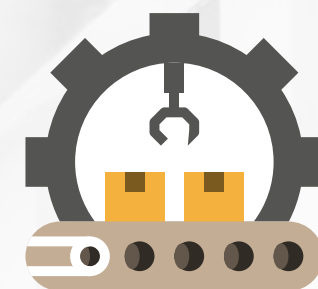
Planning in the acquisition of plant equipment

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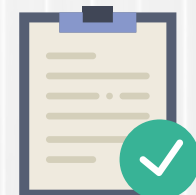
Comprehensive Analysis and Management

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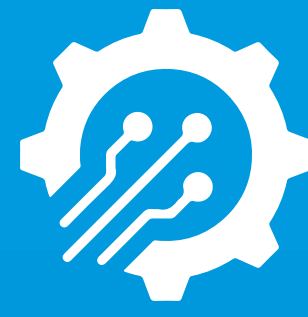
We offer comprehensive solutions, providing consultancy and training to optimize production. We design plans according to your needs.

With more than 30 years of experience behind our back, we specialize in the management of plant projects, the adaptation of different layouts, productivity improvements in GxP environments, equipment/process integration and the development of new products.

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
Process Technology

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We provide solutions for the processing of solid, semi-solid and liquid products. We specialize in the integral design and implementation of the different ways of obtaining, while applying the highest quality standards in the fulfilment of each requirement.

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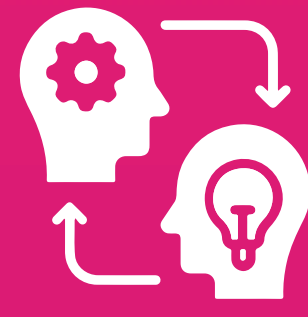Technology Transfer

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We become your strategic ally offering quality assurance in processes and products. We provide everything you need to update technological systems, guiding you through each stage in order to improve your productivity, reduce costs and facilitate an adequate scaling-up, replicating the results obtained in our laboratory to your industrial plant.

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Validations and Qualifications

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Our staff is highly trained in the preparation of supporting documentation related to the calibration, qualification and validation of production equipment. We have the necessary instruments to carry out the corresponding qualification protocols of both machines and their processes. FAT, SAT, IQ, OQ, PQ, SV (Software Validation) and CV (Cleaning Validation).

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After-sales Service

Servicio Post venta

Our After Sales department is formed by a dedicated and skilled team of mechanical, electronic, programming and administrative technicians who are dedicated to providing a fast and accurate response to our customers around the world.

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Analysis and Integral Management

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We understand the needs of our clients, devising a plan that allows us to monitor and identify their production capacities in order to minimize their costs, improve the quality control of their products, optimize efficiency (OEE) and increase productivity in their manufacturing plant.

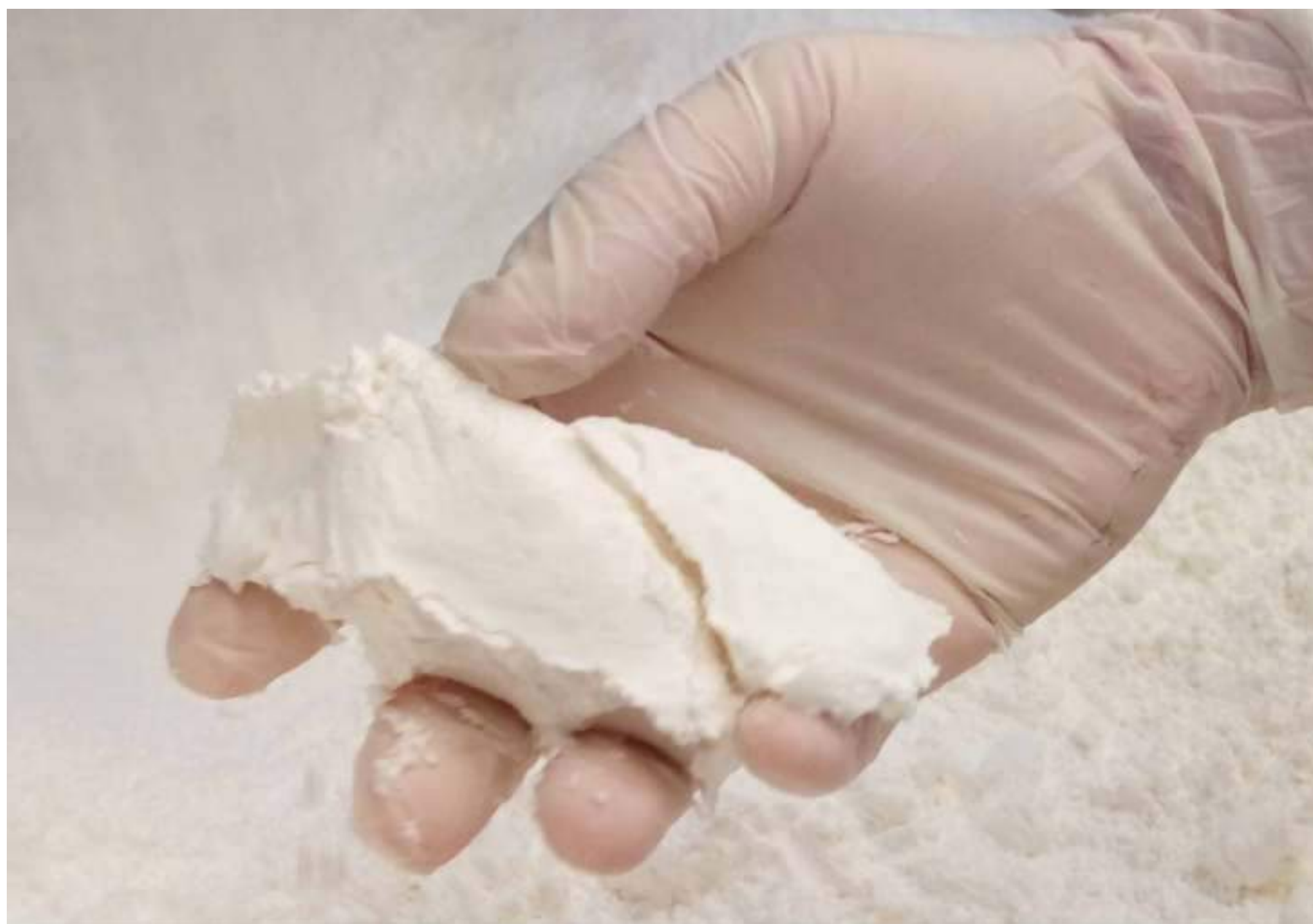
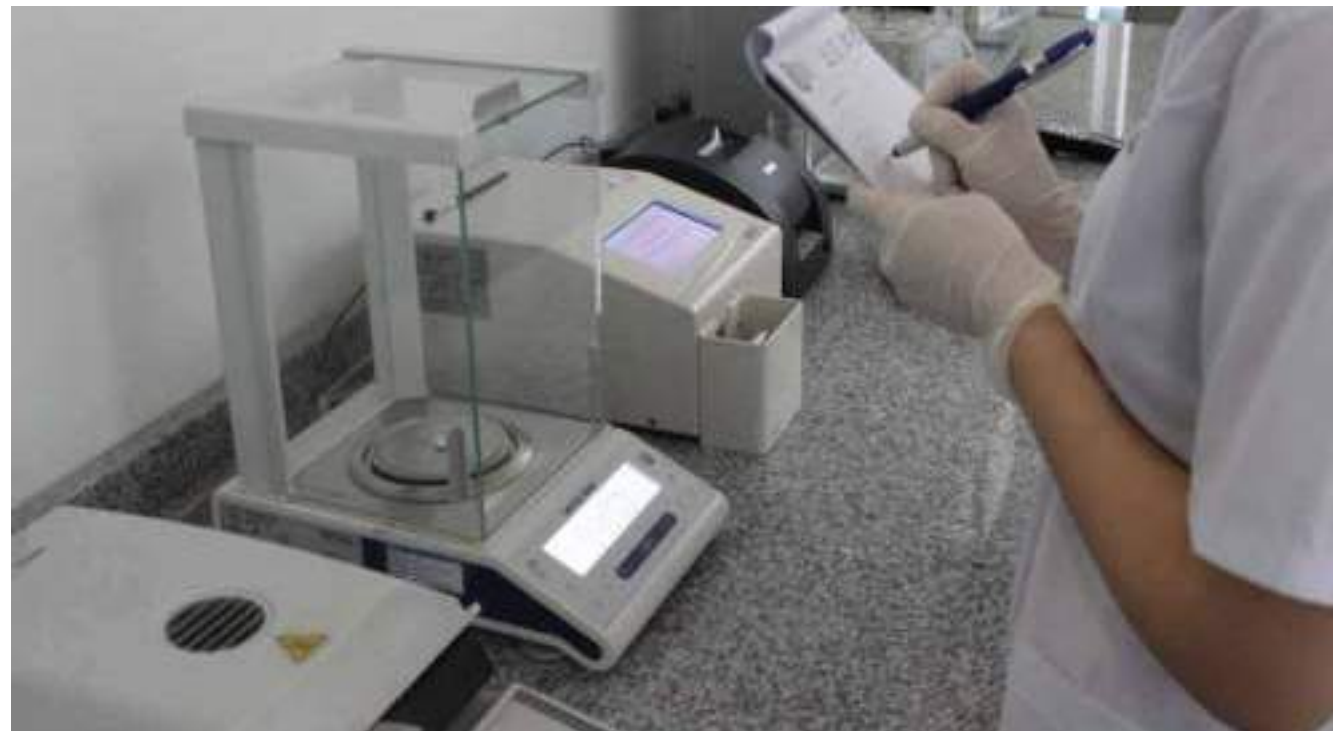
We train your human capital throughout the entire production process, detecting opportunities for improvement in each of these instances.

In this way, we automate each of these stages achieving greater agility in the supply chain, a substantial reduction in costs, manufacturing times and efficient management of your resources to maximize profitability.

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We offer equipment for solid, semi-solid and liquid lines. We have machinery for R&D accompanied by our production equipment, in small and large volumes, which allows geometric, kinematic and dynamic scaling.

We have the development to automatically wash and dry equipment, achieving a reduction in water consumption through safe and repetitive processes with the purpose of obtaining a cleaning validation.

To complete our extensive line of assistance, we make available the experience and know-how of our chemists and pharmacists, supporting our clients in the training and adaptation stage of their production processes.

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

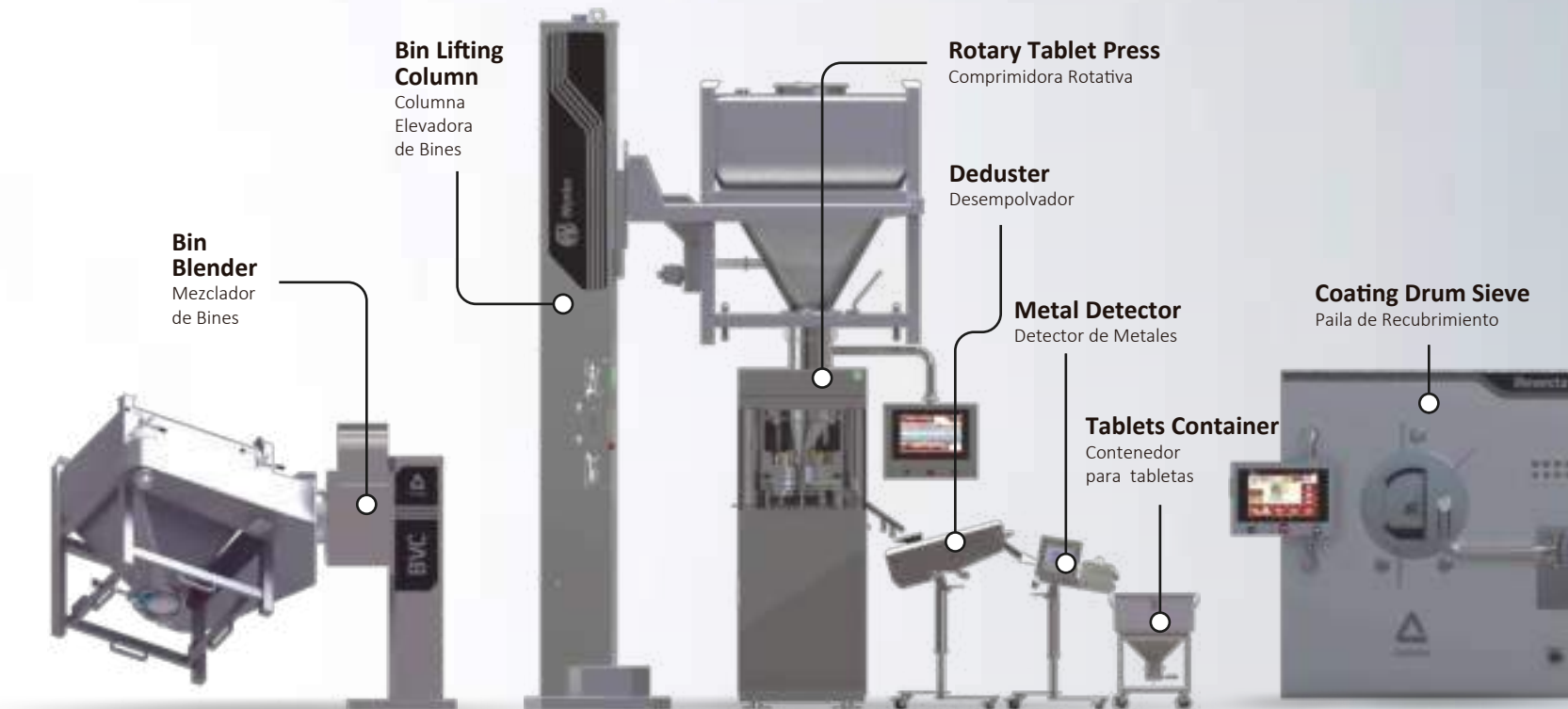
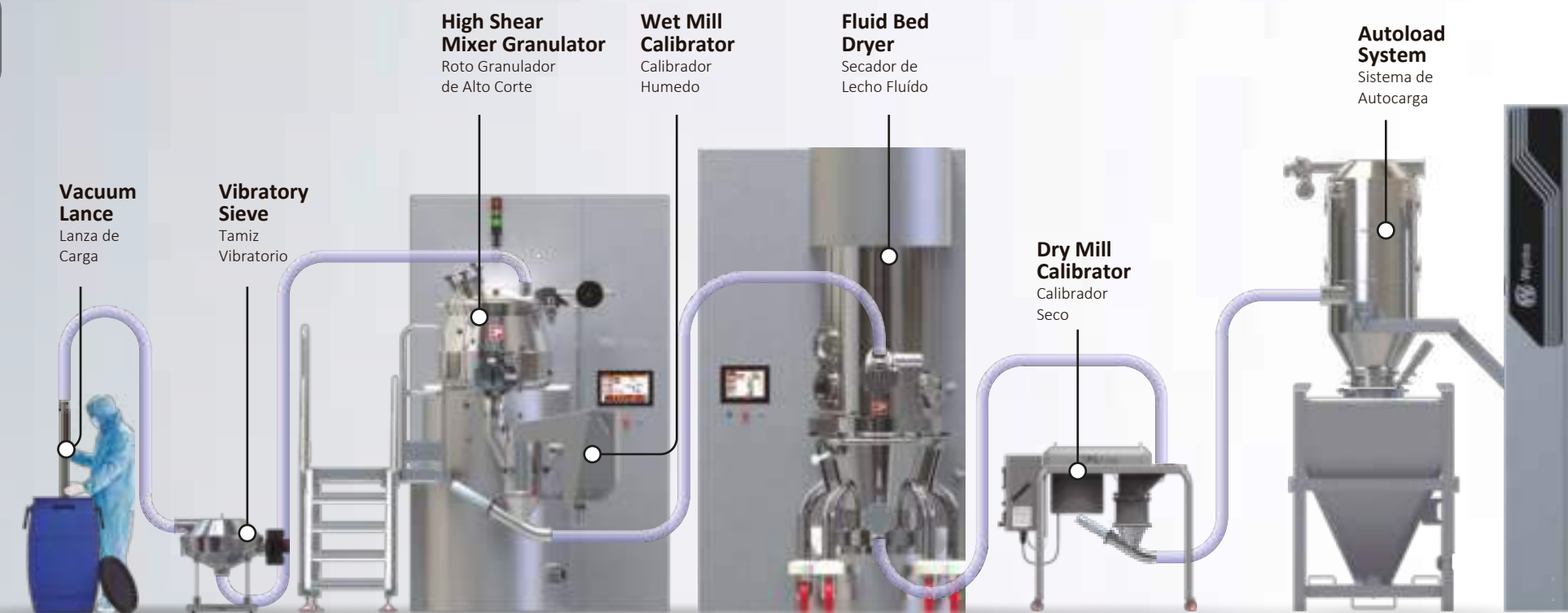
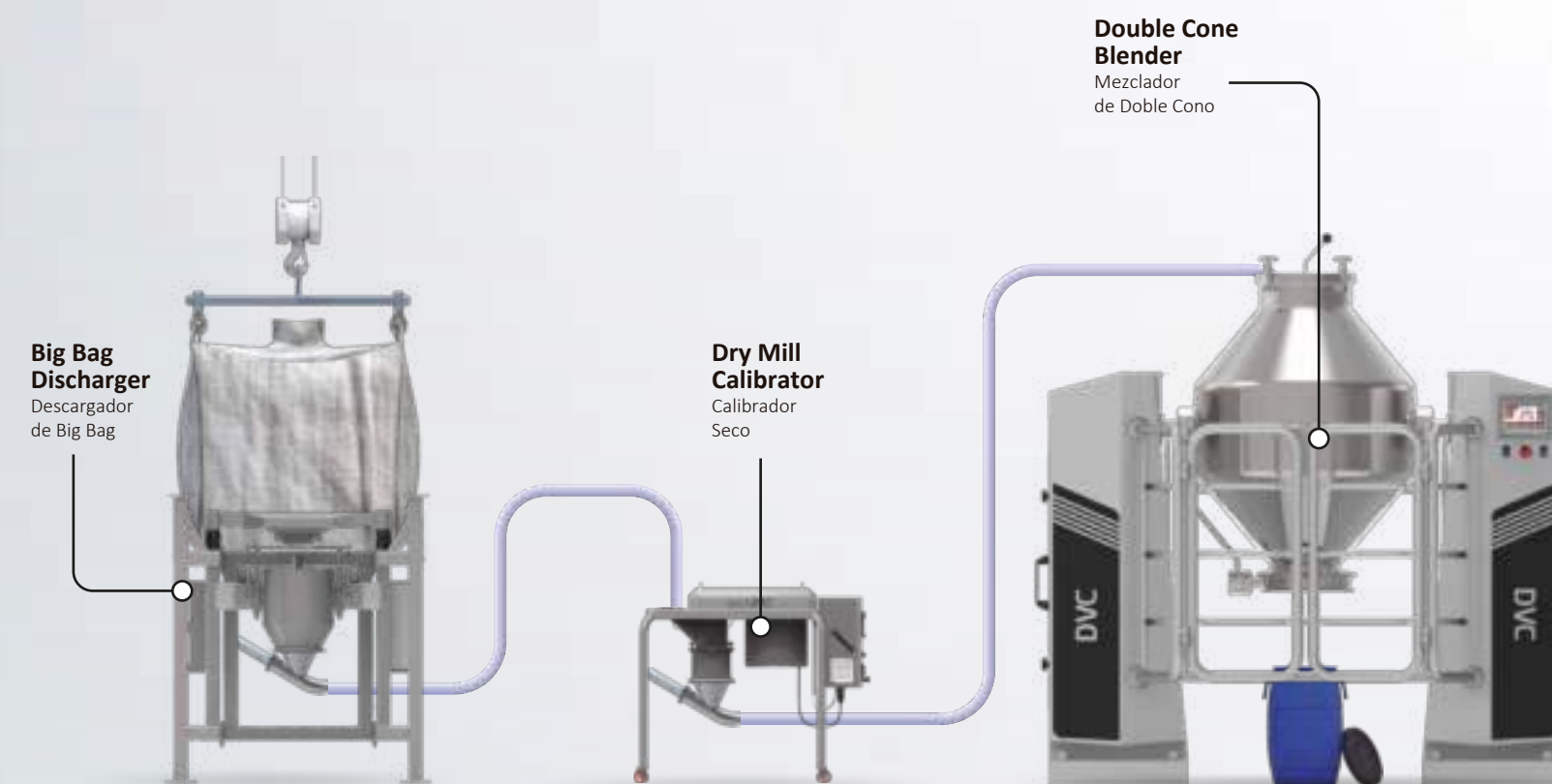
Compresión Directa

Dry Granulation

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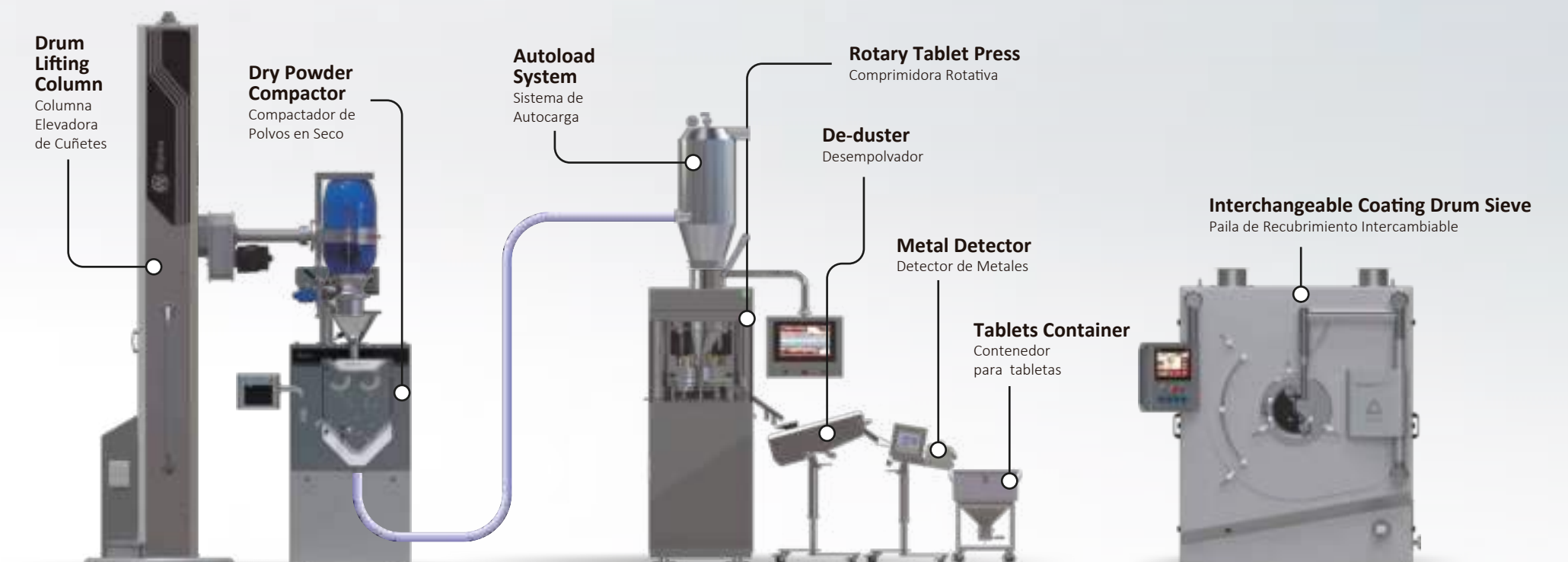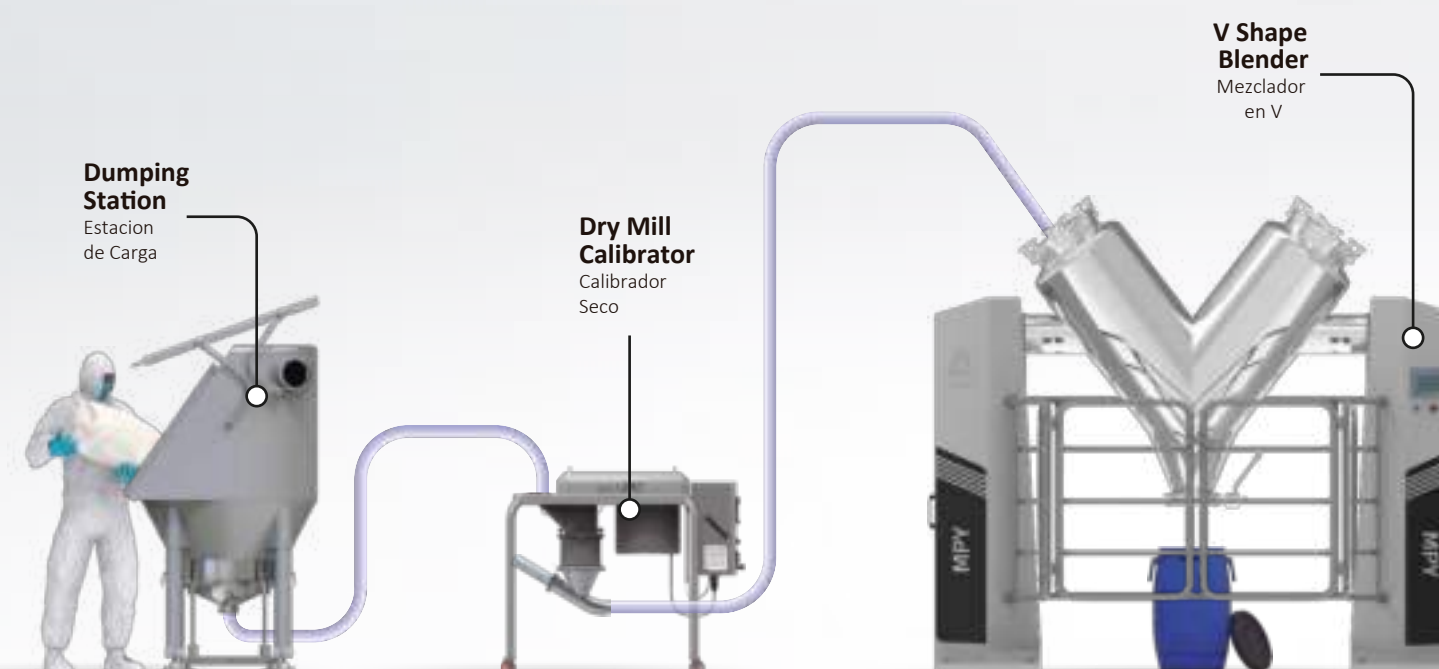MIC high shear mixer granulators are equipped and instrumented to control critical variables that regulate granulation. The wide range of mixing speed, a sensitive flow control of the liquid to be incorporated, different drop dispersing nozzles and the process cut by time or by torque, allow a precise adjustment of the granulation process.

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Development model

ESSICCA fluid bed dryer introduces two very important practical concepts:

- **Optimization of the drying process: Defines the optimum air flow for a granulate based on the determination of the VMF (minimum fluidization speed).**
- **Automation of the drying process: It makes the drying process independent of the operator through recipes. It allows estimating the end point of drying by monitoring the temperature of the product, after the drying phase at a constant speed.**

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Development model

The CPS and CPM conical calibrators are impact and shear mills, which allow precise control of the variables of the solid particle calibration process (size reduction) both in raw materials and in dry and wet granules.

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Sizes | Medidas

4x4mm
6x6 mm
8x8 mm
10x10mm
12x12 mm
15x15 mm

The BVC bin blender has been conceived taking into account the three main blending variables:

- A) The geometry of the container.
- B) The filling level or load of the equipment.
- C) The blending speed.

Powder mixes with "free flow" characteristics can be blended for a given time regardless of the chosen speed (RPM), while those of the "cohesive" type must be blended at high speeds to facilitate the shearing and mixing of the components.

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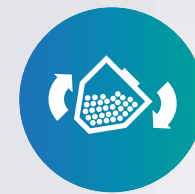
2000 liters

Double Cone Blender

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The double cone blender DVC Pharma has been conceived taking into account the three main blending variables:

- A) The geometry of the container.
- B) The filling level or load of the equipment.
- C) The blending speed.

Powder mixes with "free flow" characteristics can be blended for a given time regardless of the chosen speed (RPM), while those of the "cohesive" type must be blended at high speeds to facilitate the shearing and mixing of the components.

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Development model

V-Shape Blender

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The V-shape blender MPY Pharma has been conceived taking into account the three main blending variables:

- A) The geometry of the container.
- B) The filling level or load of the equipment.
- C) The blending speed.

Powder mixes with "free flow" characteristics can be blended for a given time regardless of the chosen speed (RPM), while those of the "cohesive" type must be blended at high speeds to facilitate the shearing and mixing of the components.

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Lifting Columns

Bin, drum and autoload lifting columns

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The design of our columns can be used with other pharmaceutical equipment such as blenders, coaters, tablet presses, capsule filling machines, etc.

They are available in different options with unique and modern technologies to optimize operations in the pharmaceutical industry.

Depending on the customer's requirement, the lifting column will be designed to transfer or unload powders, granules, tablets or other products through the use of drums, bins or autoload systems.

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The Kompressor tablet press, developed under an innovative concept, allows format change in an agile and safe way. Its structural robustness makes it possible to work in difficult compression condition during prolonged periods.

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Kompressor 200

Kompressor 300

Kompactor is a powder compactor designed for dry granulation, especially for pilot batches and production, which meets the highest requirements and standards of the pharmaceutical industry.

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~200 kg / h

Development model

DS6000 is a tablets dust remover equipped with a powder aspiration system and helical advance screw. Its simplicity and compact design make it an affordable plug-and-play product.

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DT8000 is a metal detector and separator with maximum sensitivity for detecting metals ferrous type and nonferrous, as well as stainless steel.

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Revesta is a pail designed to perform coating on a galenic scale or production batches of pellets or tablets, and can be used in both aqueous and organic coatings, ensuring product quality and reproducibility.

It is equipped with different levels of alarms, this completes a configuration that provides safety to the operator in a process as critical as coating.

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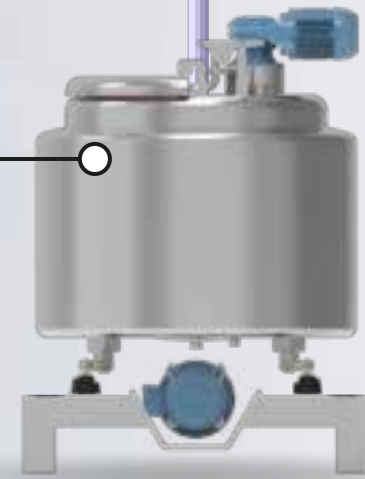
Development model


Semi-solids Production Line with Packaging Storage Tank

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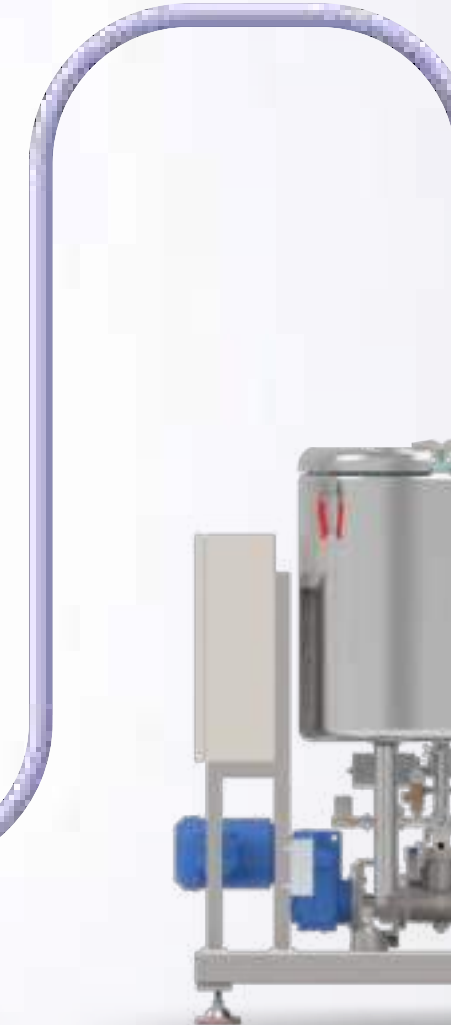
Storage Tank
Tanque para Almacenamiento

To Filling Machine
A Envasadora


Semi-solids Production Line with Mobile Reactor

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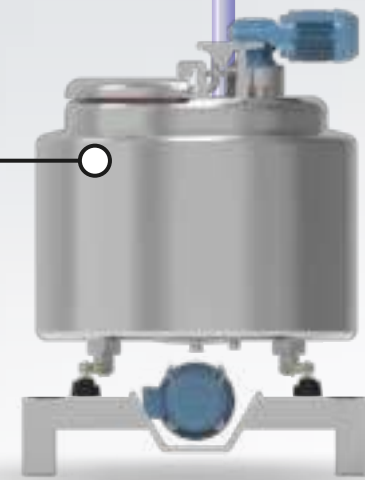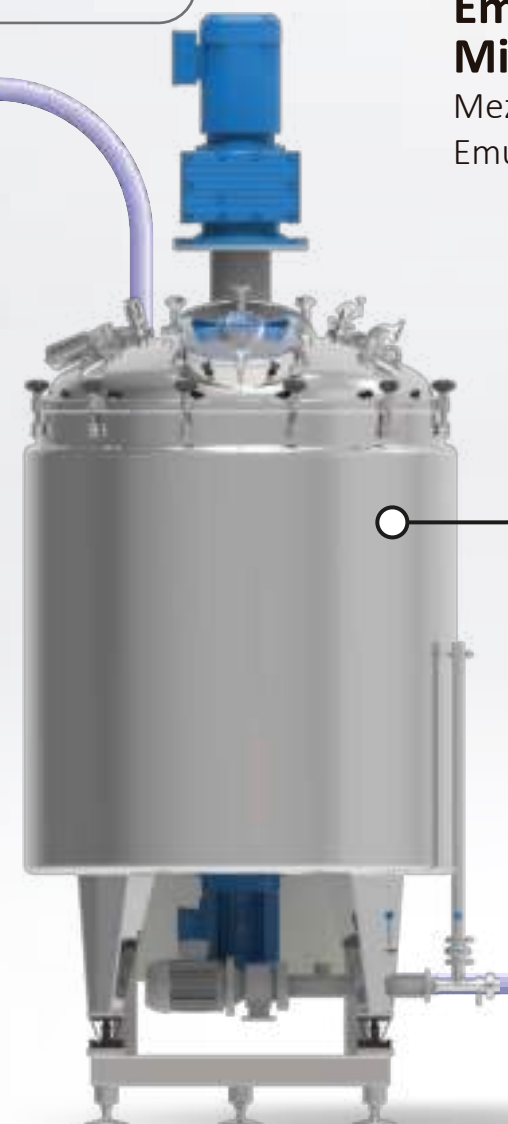
Transfer Reactor
Reactor para Trasvasado

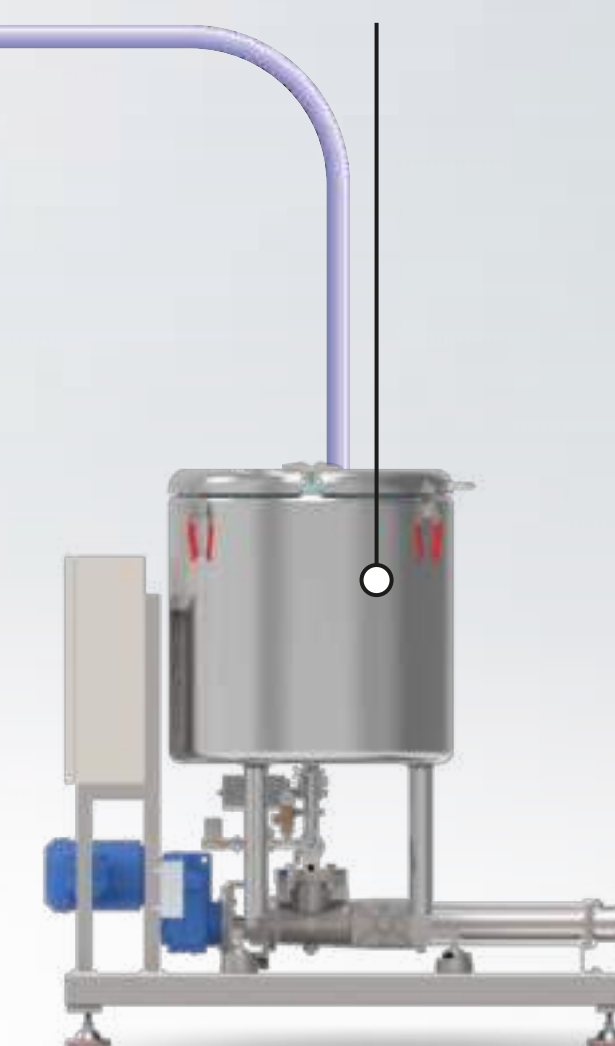Storage Tank
Tanque para Almacenamiento

To Filling Machine
A Envasadora

turbomixer designed for the manufacture of creams and gels, achieving very short mixing and emulsifying times. In its development version, it allows the production of small reproducible batches on a productive scale, fundamentally due to its geometric, kinematic and dynamic similarity.

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Development model

Designed for the production of creams and gels with very short mixing and emulsifying times, this single control reactor is used for the production of medium and high viscosity products.

Among its main features, it has anchor type agitation with scrapers together with the possibility of incorporating an emulsifier.

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
Liquid Production Line

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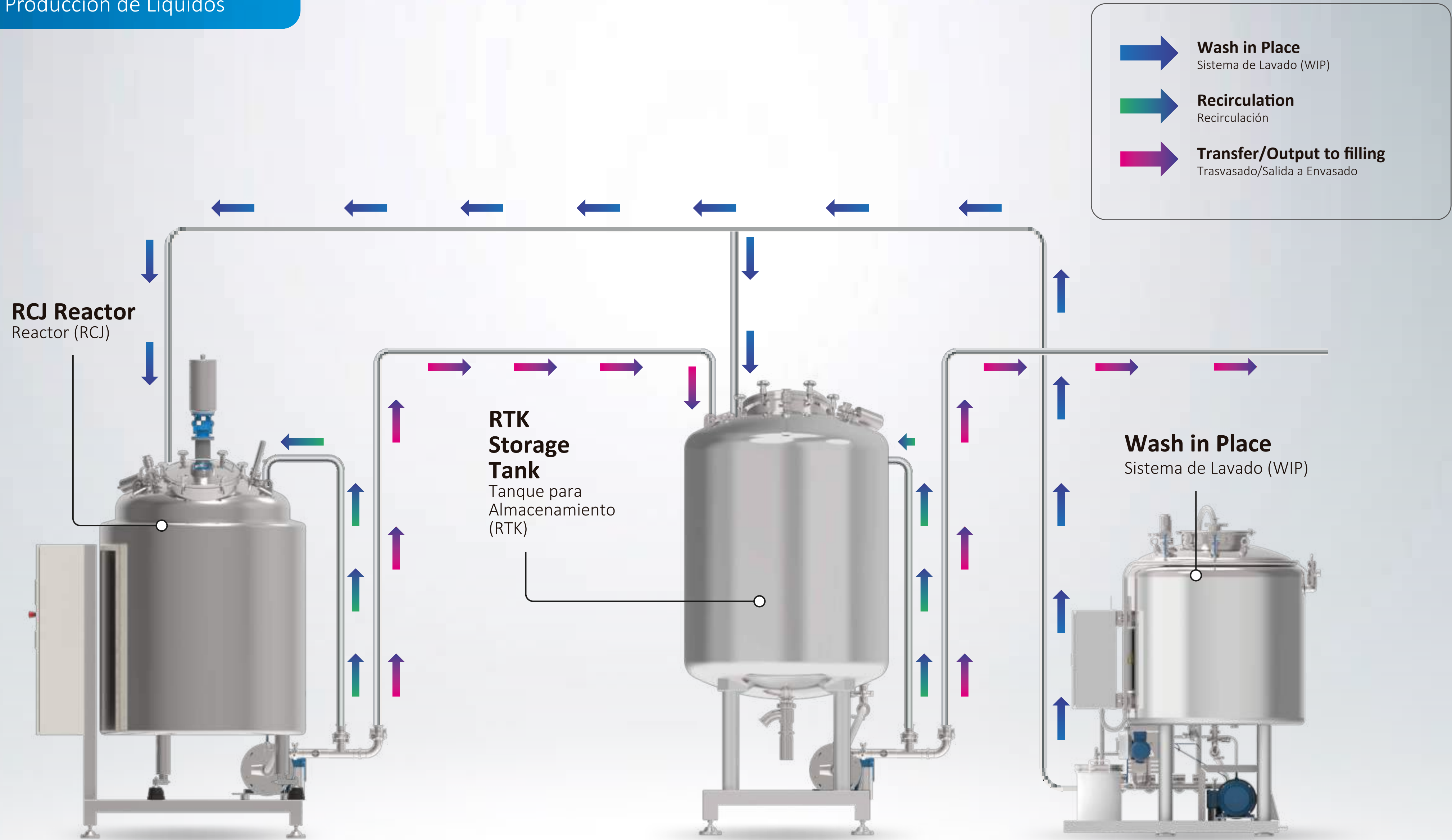Reactors

Different reactors type for multiple uses

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Our line of reactors includes sterile products, syrups, suspensions and biotechnology, among others. They are built in stainless steel with AISI 316L quality under GMP guidelines and have a connection for cleaning that allows the automatic washing and sanitizing process.

According to the requirements, they admit heating and cooling; with waterproof insulation. They have a design with sanitary loading and unloading valves as well as a control system according to the specific requirement of the client.

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Binders, Emulsions, Steriles, Transport, Storage and other uses

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Adjusted to pharmaceutical standards, the wash in place (WIP) system is an autonomous equipment that has the possibility of washing solid, semi-solid and liquid process lines.

Designed under GMP guidelines, among its salient features it allows the possibility of editing automatic cleaning processes, a simple implementation in plant reducing production downtime, saving water and associated chemical agents and a simple integration to centralised process data acquisition systems.

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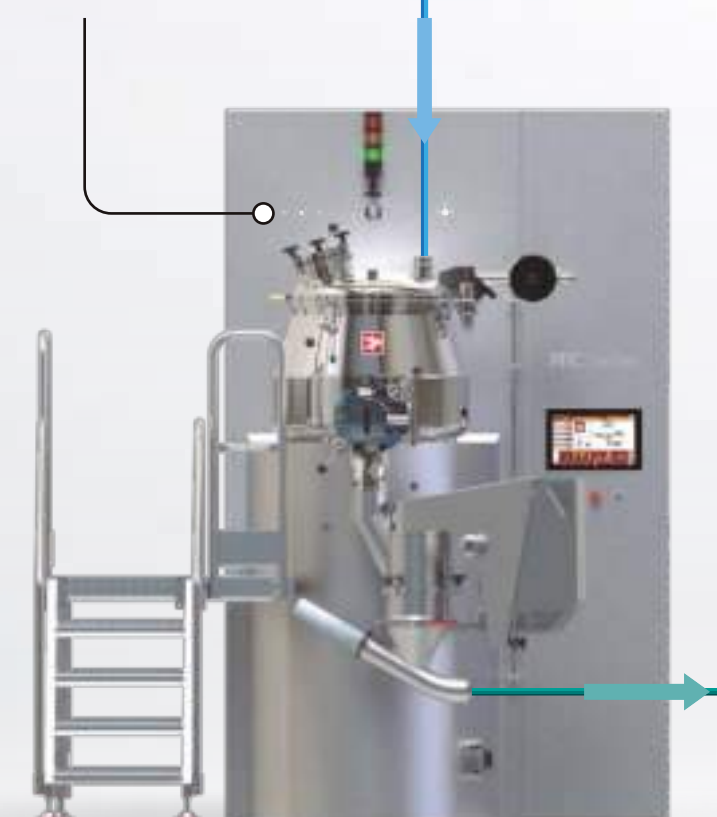
Washing of Equipment Lavado de Equipamiento

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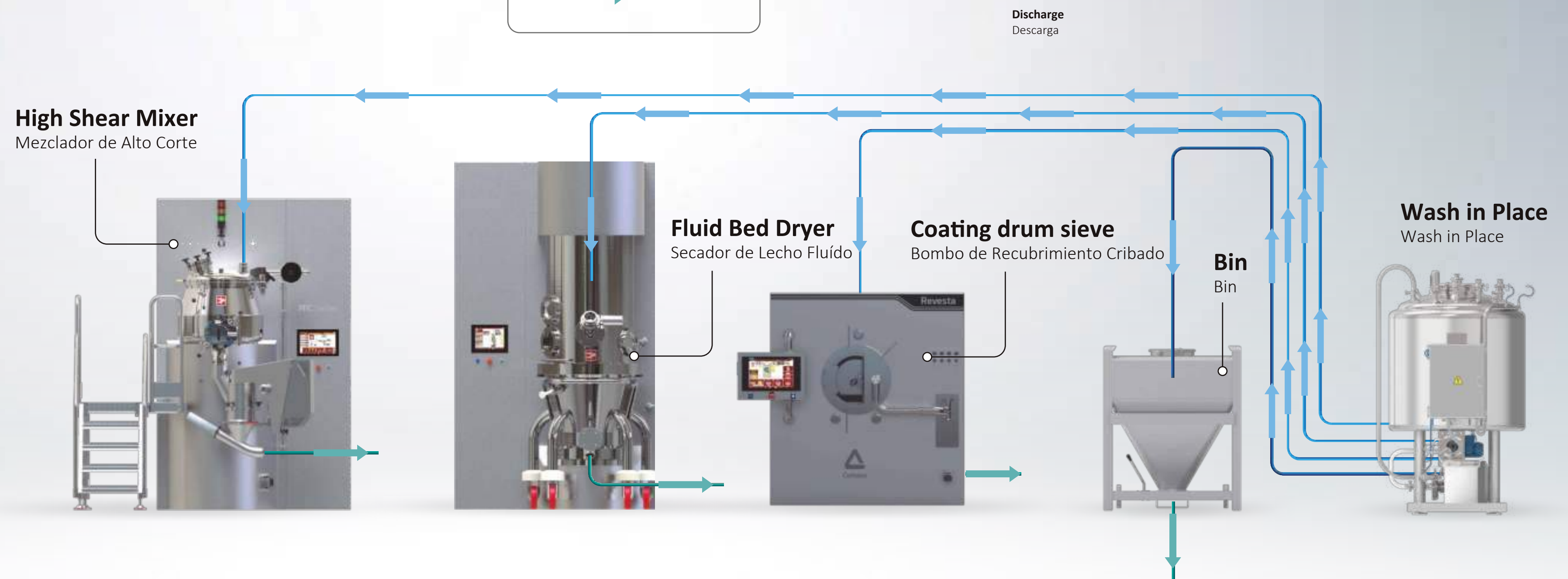
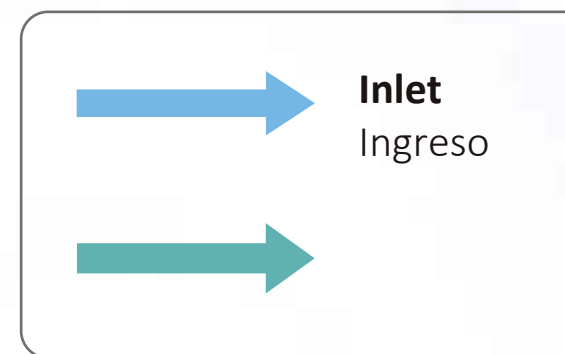
Bin Bin

Wash in Place Wash in Place


This system is based on a sanitary air treatment adapted to the solids line by means of thermal transfer by hot air, thus optimising labour times in this operation.

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
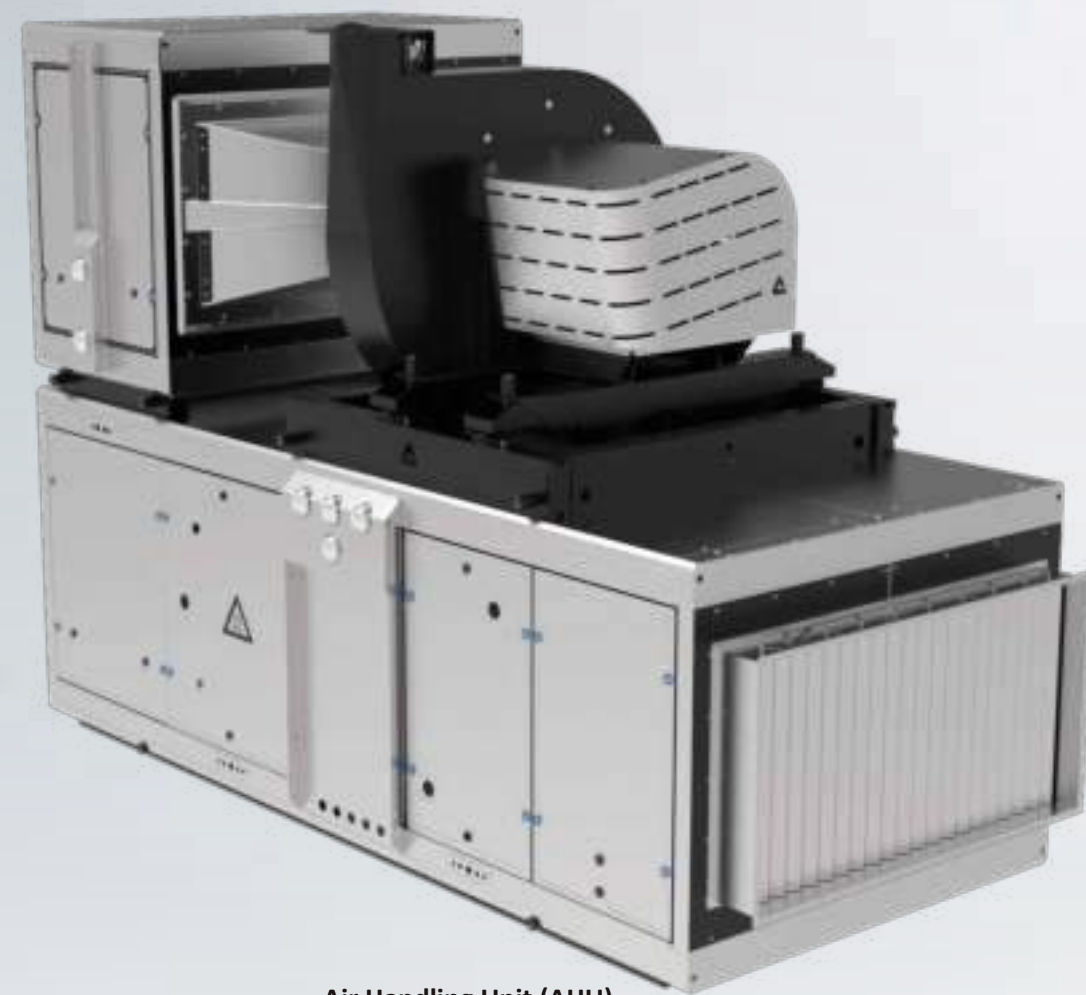
Drying of Equipment
Secado de Equipamiento

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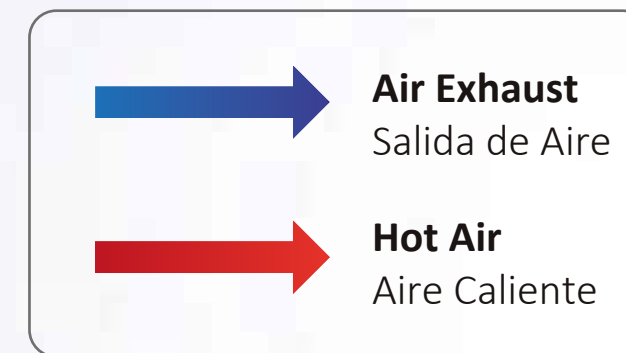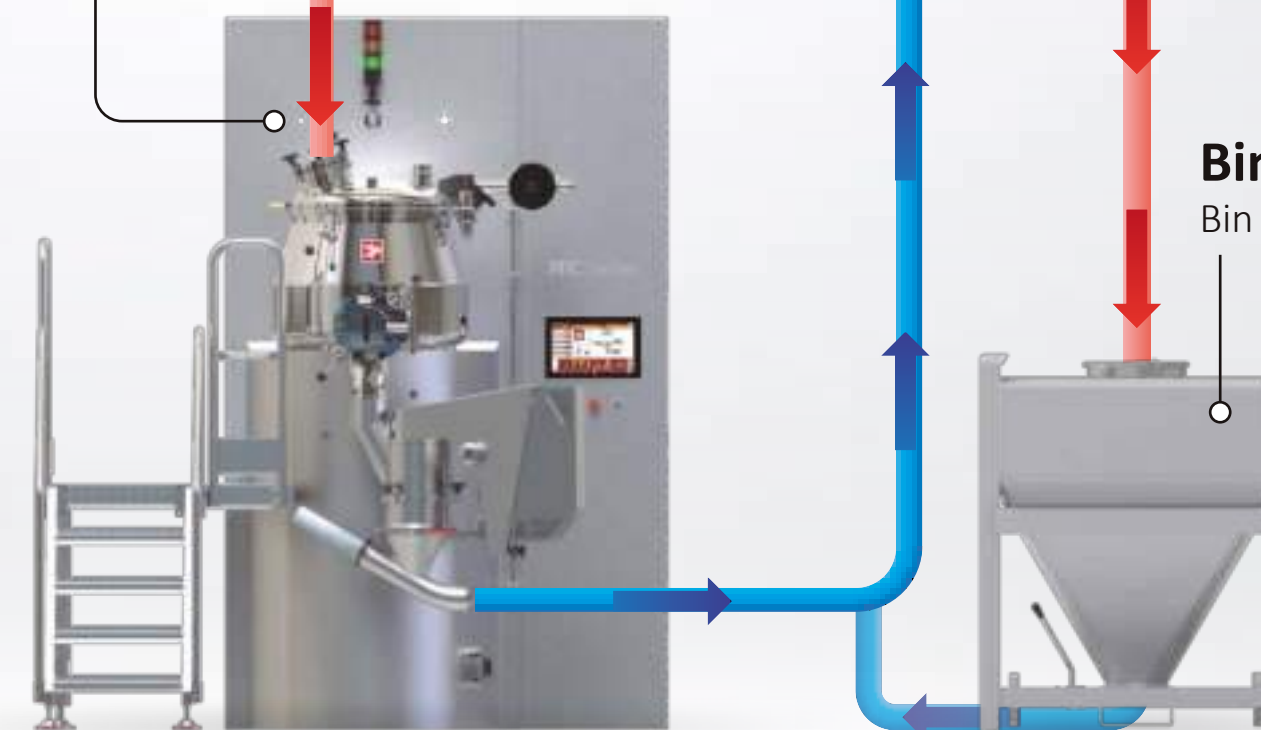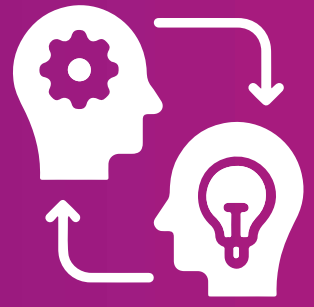
Technology Transfer

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We have a laboratory for new product development and customer support. We have the latest technology in terms of equipment and analytical instruments that allow us to carry out process controls during each stage, from formulation to the delivery of documentation for registration.

This is complemented by our outstanding team of professionals who have extensive experience in the industry, with a large number of products developed in various solid, semi-solid and liquid dosage forms.

In this way, our service generates an extensive cost reduction and avoids the so-called "trial and error", avoiding incurring additional expenses and delays to start production with the new technology without setbacks.

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Validations and Qualifications

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Good manufacturing practices and international quality standards describe the design, manufacture, installation and operation of equipment. This implies that they must be routinely calibrated, inspected or verified according to a pre-written programme designed to ensure proper performance and compliance with GMP regulations.

The documentation represents the intellectual capital that the company must preserve in order to standardise and improve processes, carry out correct organisational planning and effective control of the activities and accredit the material required at the time of an audit.

Our team of professionals is in charge of the supporting documentation, related to the calibration, qualification and validation of production equipment, in order to prepare and execute the DQ, FAT, SAT, IQ, OQ, PQ protocols.

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After-Sale Services

Servicio Post Venta

When you buy our equipment, you choose much more than a machine. You opt for a work philosophy that respects principles and values, a relationship that begins by placing your trust in us.

Based on the philosophy of continuous improvement, we are present on all five continents with our representatives and we invest in their training and development, offering an efficient response that allows us to provide the quality of service that our customers deserve.

We maintain an exclusive and individual treatment through the provision of original spare parts, accompanied by a review and diagnosis service that preserves their optimal operating and safety conditions. In this way, we prolong their useful life, backed by an extended three-year warranty and a hardware upgrade to update equipment that is more than 10 years old.

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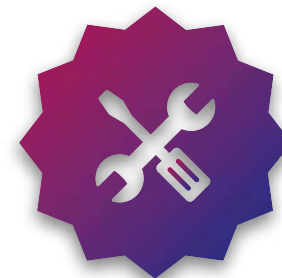
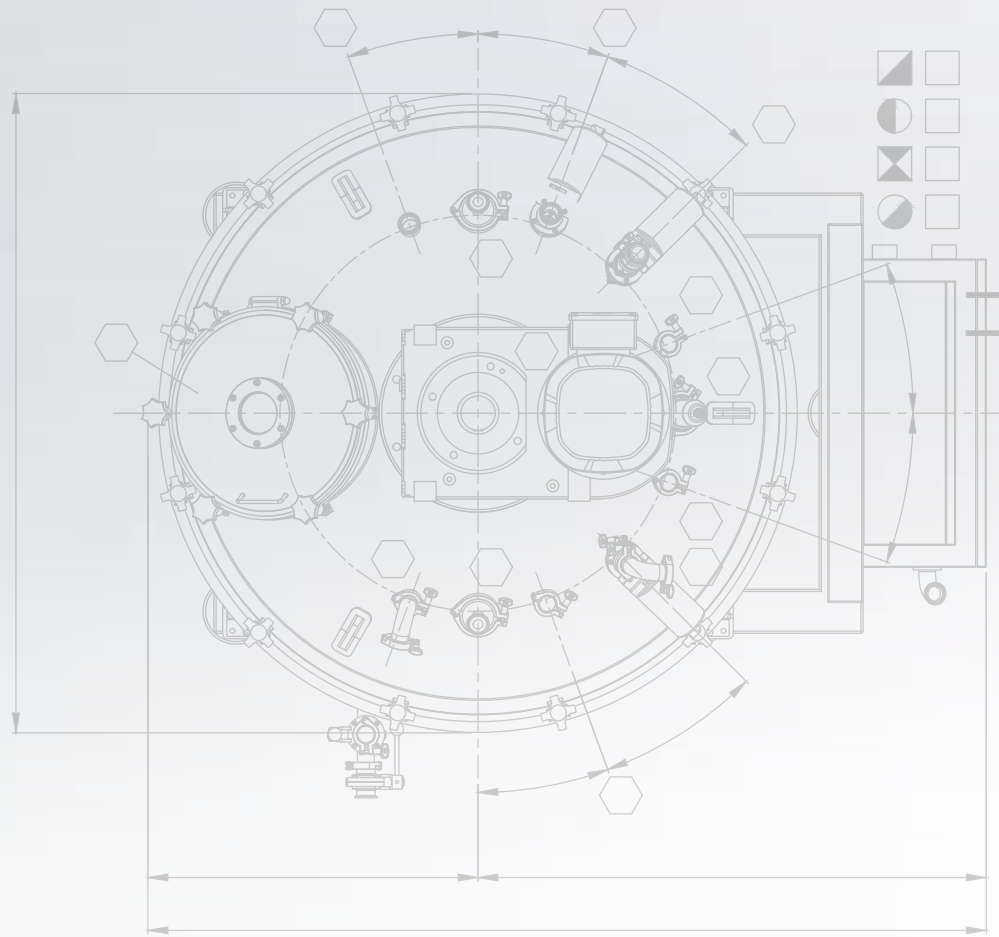
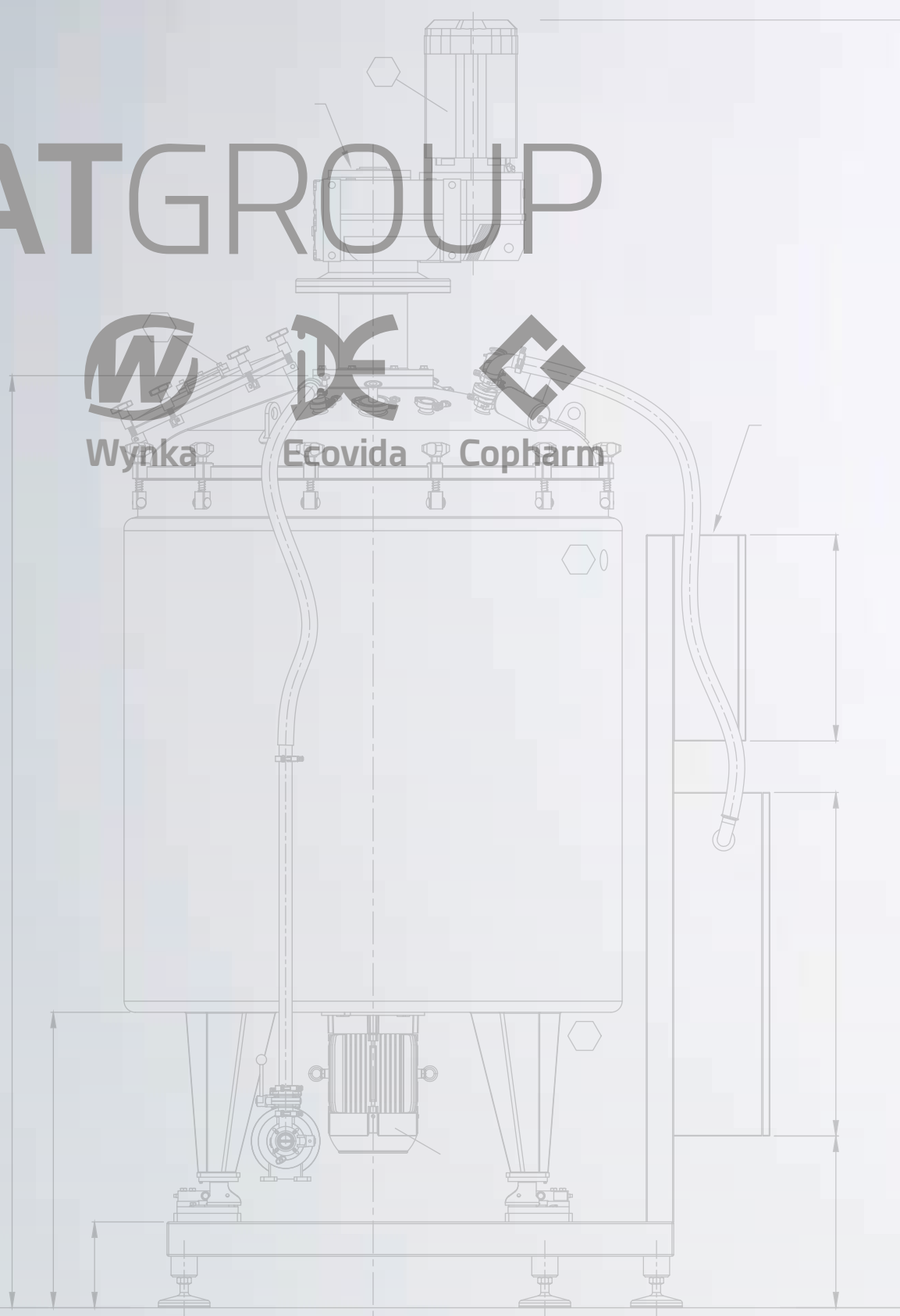
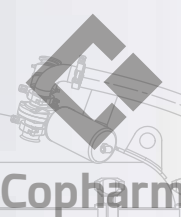
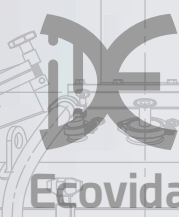
 **Responsibility**
Responsabilidad

PATGROUP

Built in compliance with GMP guidelines

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Software in compliance with 21 CFR Part 11 regulations

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Equipment in compliance with ATEX and EX-PROOF regulations

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Batch Record in compliance with 21 CFR Part 11 regulations

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